

# NanoQAM

*Centre de Recherche sur les Nanomatériaux et l'Énergie*

## ***QUICK START***

### **Hot Cathode Ionization Vacuum Gauge**

*By The Hornet InstruTech*

## How to start the high vacuum gauge



- 1- Press menu button
- 2- Chose setup unit + press enter
- 3- Chose FP operate + press enter
- 4- Chose unit on/off + press enter
- 5- Using the arrows switch from off to on + press enter

If you have any question, please contact Galyna Shul ([shul.galyna@uqam.ca](mailto:shul.galyna@uqam.ca), 514 250 5306).