

	SERVICE REFERENCE		
	Product(s):	Advantage & QNX	Prepared By:
	Subject:	<i>How to Perform Reset Saved Parameters from the QNX Touchscreen.</i>	

1. Close Advantage
2. Right click on the **Windows Taskbar** and select **Start Task Manager**.
3. After **Task Manager** opens, enable the checkbox “**Show processes from all users**”.



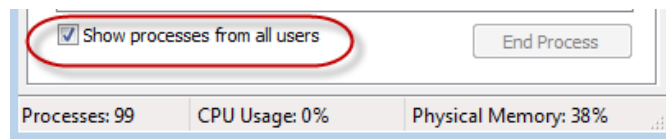

4. Now select the **Processes** tab and verify that the **QAdvX.exe** process for the instrument that requires the **Reset Parameters** is not listed as running. If the process is listed, highlight the process and then right click and select **End Process** to terminate the application.

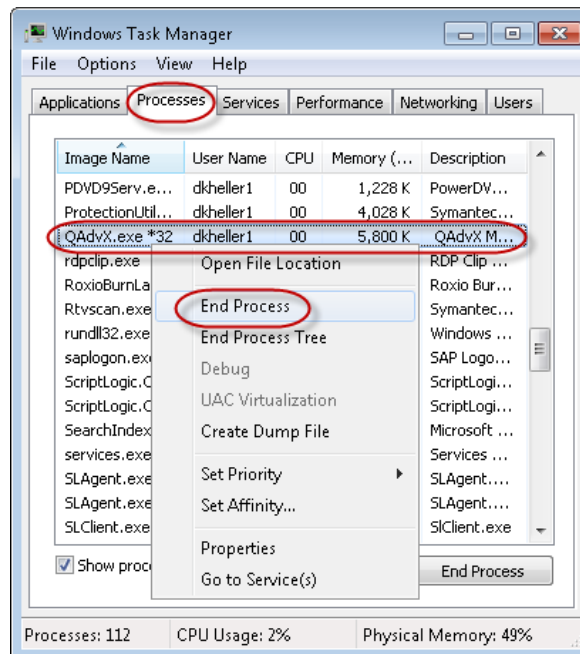


5. After the processes **QAdvX.exe** has been terminated, go to the Instrument touch screen.
6. Press the **Control** menu button.




	SERVICE REFERENCE		
	Product(s):	Advantage & QNX	Prepared By:
	Subject:	<i>How to Perform Reset Saved Parameters from the QNX Touchscreen.</i>	


7. Now click the white down triangle  until the menu item Reset Saved Parameters appears.

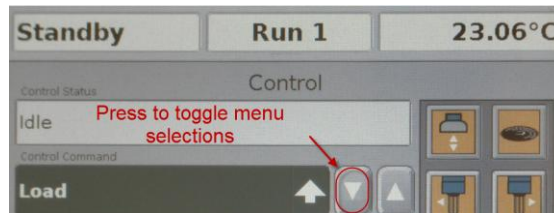


8. Now press the menu item **Reset Saved Parameters**.

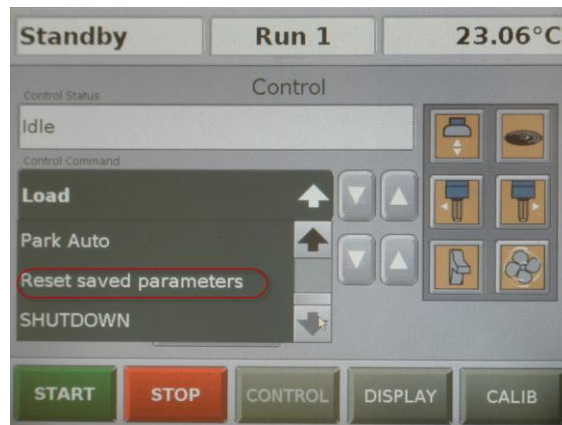


9. Now press the Apply button.

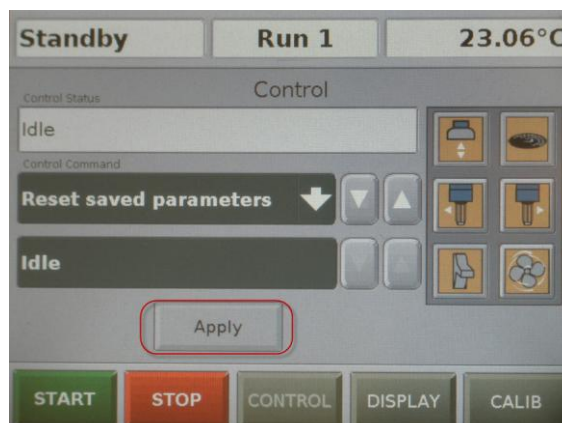



	SERVICE REFERENCE		
	Product(s):	Advantage & QNX	Prepared By:
	Subject:	<i>How to Perform Reset Saved Parameters from the QNX Touchscreen.</i>	

10. When the following message appears on the screen, click the **YES** button.

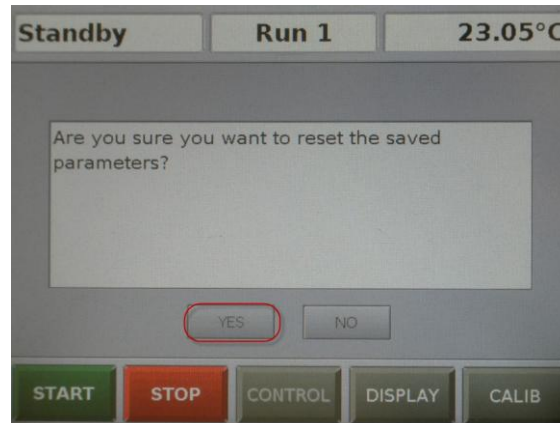


11. The following message, #735 may also appear during this process depending on the instrument type. If the message appears on the screen, click the **OK** button.

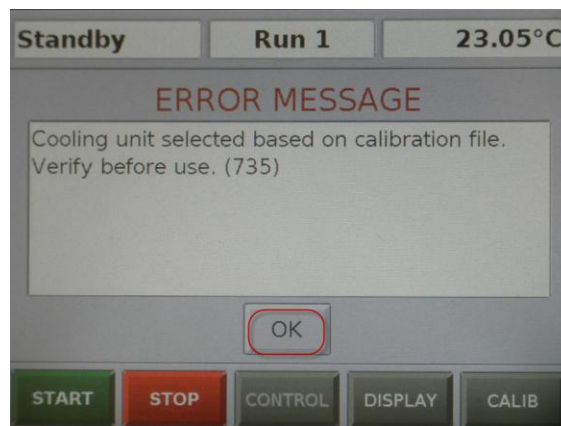

	SERVICE REFERENCE		
	Product(s):	Advantage & QNX	Prepared By:
	Subject:	<i>How to Perform Reset Saved Parameters from the QNX Touchscreen.</i>	

12. When the following message, #977 appears on the screen, indicating that the **Instruments Parameters** have been successfully reinitialized, click the **OK** button.

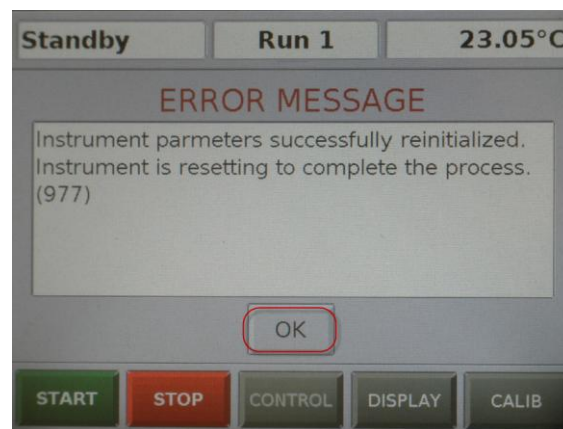

13. The instrument will now automatically reboot.

14. When the instrument has completed rebooting, the TA logo will now appear on the touch screen.

15. Now connect to the instrument by going to TA Explorer and double clicking on the icon for your instrument.